

## DAB Differentiating Solution Kit

### Contents:

#### Volume

125 ml DAB Differentiating Solution 1  
125 ml DAB Differentiating Solution 2  
250 ml Decolorizer

**Storage:** Store in dark at room temperature.

### Procedure:

1. To prepare working solution, mix DAB Differentiating Solution 1, DAB Differentiating Solution 2 and DI water in the following ratio. Mix well (Note: solution will be purple).  

98% DI water  
1% DAB Away 1  
1% DAB Away 2
2. Apply to stained tissue section.
3. Allow to incubate 1-10 minutes.
4. Rinse with water.
5. Apply Decolorizer to surfaces for 5-10 seconds.

**Precautions:**

- \* All three components are acidic.
- \* Avoid contact with skin or eyes.
- \* Wear protective gloves and eye covers.
- \* In case of contact, wash area with copious amounts of water.