

# Normal Antibody Diluent (Phosphate Buffered)

**Description:**

Provided as a ready-to-use (Phosphate Buffered) diluent for diluting primary antibodies and for use as a negative control. Both monoclonal and polyclonal antibodies are stabilized for long-term storage thereby reducing the number of titrations required from concentrated form. This product contains a highly purified carrier protein that minimizes proteolytic enzyme degradation, a wetting agent to reduce surface tension and an anti-microbial agent to prevent microbial growth. Not for antibodies that have been conjugated with peroxidase. In most cases, antibodies diluted in this reagent can be stored for 24 months at 2-8°C. Diluent is subjected to 0.2 micron filtration.

pH 7.4

**Availability/Contents:**

<u>Item #</u>	<u>Volume</u>
ABB125	125 ml
ABB500	500 ml
ABB999	1000 ml

**Uses/Limitations:**

Not to be taken internally.  
For In-Vitro Diagnostic use.  
Histological applications.  
Do not use if reagents become cloudy.  
Do not use past expiration date.  
Use caution when handling reagents.  
Non-Sterile.



Ordering Information and Current Pricing at [www.scytek.com](http://www.scytek.com)

**Storage:**

Store at 2-8°C.


**Precautions:**

Avoid contact with skin and eyes.  
Harmful if swallowed.  
Follow all Federal, State, and local regulations regarding disposal.

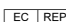
**Procedure:**

1. Pour appropriate volume of Primary Antibody Diluent into clean mixing bottle.
2. Dilute primary antibody directly in Primary Antibody Diluent (Phosphate, Green).
3. Use as directed by antibody manufacturer.
4. Store diluted antibody at 2-8°C.

Storage: 2° C  8° C

 ScyTek Laboratories, Inc.  
205 South 600 West  
Logan, UT 84321  
U.S.A.


 

  
Emergo Europe  
Prinsessegracht 20  
2514 AP The Hague, The Netherlands

### References:

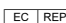
1. Contreras GA, Thelen K, Ayala-Lopez N, Watts SW. The distribution and adipogenic potential of perivascular adipose tissue adipocyte progenitors is dependent on sexual dimorphism and vessel location. *Physiological reports*. 2016 Oct 1;4(19):e12993.
2. M. A. Bates, C. Brandenberger, I. Langohr, K. Kumagai, J. R. Harkema, A. Holian, and J. J. Pestka, "Silica Triggers Inflammation and Ectopic Lymphoid Neogenesis in the Lungs in Parallel with Accelerated Onset of Systemic Autoimmunity and Glomerulonephritis in the Lupus-Prone NZBWF1 Mouse," *PLoS ONE*, vol. 10, no. 5, p. e0125481, May 2015.
3. Kopec AK, Joshi N, Towery KL, Kassel KM, Sullivan BP, Flick MJ, Luyendyk JP. Thrombin Inhibition with Dabigatran Protects against High-Fat Diet–Induced Fatty Liver Disease in Mice. *Journal of Pharmacology and Experimental Therapeutics*. 2014 Nov 1;351(2):288-97.
4. Ono H, Basson MD, Ito H. PTK6 promotes cancer migration and invasion in pancreatic cancer cells dependent on ERK signaling. *PloS one*. 2014 May 1;9(5):e96060.
5. Wehrwein EA, Novotny M, Swain GM, Parker LM, Esfahanian M, Spitsbergen JM, Habecker BA, Kreulen DL. Regional changes in cardiac and stellate ganglion norepinephrine transporter in DOCA–salt hypertension. *Autonomic Neuroscience*. 2013 Dec 31;179(1):99-107.
6. Karmaus PW, Chen W, Crawford RB, Harkema JR, Kaplan BL, Kaminski NE. Deletion of cannabinoid receptors 1 and 2 exacerbates APC function to increase inflammation and cellular immunity during influenza infection. *Journal of leukocyte biology*. 2011 Nov 1;90(5):983-95.
7. White MG, Wang Y, Akay C, Lindl KA, Kolson DL, Jordan-Sciutto KL. Parallel high throughput neuronal toxicity assays demonstrate uncoupling between loss of mitochondrial membrane potential and neuronal damage in a model of HIV-induced neurodegeneration. *Neuroscience research*. 2011 Jun 30;70(2):220-9.
8. Swanlund, J.M., Kregel, K.C., Oberley, T.D. Investigating Autophagy, Quantitative morphometric analysis using electron microscopy. *Autophagy*, February 2010; 6(2): pages 270-277.
9. Mackenzie, C.D., Huntington, M.K., Wanji, S., Lovato, R.V., Eversole, R.R., Geary, T.G. The Association of Adult *Onchocerca volvulus* with Lymphatic Vessels. *Journal of Parasitology*. February 2010, Volume 96, Number 1, pages 219-221.
10. Oberley, T.D., Swanlund, J.M., Zhang, H.J., Kregel, K.C. Aging Results in Increased Autophagy of Mitochondria and Protein Nitration in Rat Hepatocytes Following Heat Stress. *Journal of Histochemistry and Cytochemistry*. June 2008, 56(6): pages 615-627.
11. van Maldegem, F., de Wit, M., Morsink, F., Musler, A., Weegenaar, J., van Noesel, C.J. Effects of Processing Delay, Formalin Fixation, and Immunohistochemistry on RNA Recovery From Formalin-fixed Paraffin-embedded Tissue Sections. *Diagnostic Molecular Pathology*, March 2008, Volume 17, Issue 1, pages 51-58.
12. Elstrodt, F., Hollestelle, A., Nagel, J.H.A., Gorin, M., Wasielewski, M., van den Ouweland, A., Merajver, S.D., Ethier, S.P., Schutte, M. BRCA1 Mutation Analysis of 41 Human Breast Cancer Cell Lines Reveals Three New Deleterious Mutants. *Cancer Research*, January 2006, Volume 66, Issue 1, pages 41-42.
13. Williams, S.M., Fitzgerald, S.D., Willie, M.R., Lee, L.F., Fadly, A.M. Tissue Tropism and Bursal Transformation Ability of Subgroup J Avian Leukosis Virus in White Leghorn Chickens. *Avian Diseases*, 2004, 48: pages 921-927.
14. Oberley, T.D., Zhong, W., Szweda, L.I., Oberley, L.W. Localization of Antioxidant Enzymes and Oxidative Damage Products in Normal and Malignant Prostate Epithelium. *The Prostate*, 2000, 44: pages 144-155.

 Storage: 2° C  8° C



ScyTek Laboratories, Inc.  
205 South 600 West  
Logan, UT 84321  
U.S.A.

  
Emergo Europe  
Prinsessegracht 20  
2514 AP The Hague, The Netherlands